
AccuPoint[®]

AccuPoint Advanced Sanitation Verification System
is accurate in detecting ATP.*


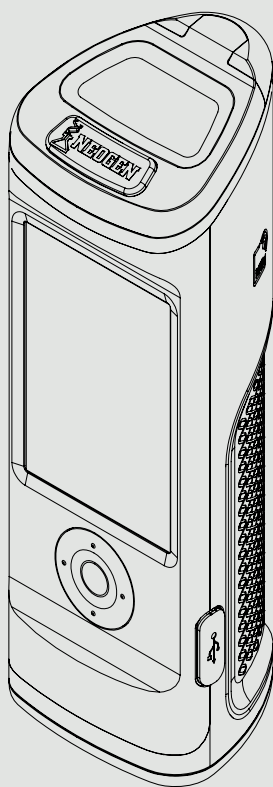
*The first sanitation verification device to be approved by the AOAC Research Institute

01. To power on the reader, press and hold the center button on your AccuPoint Phoenix. Enter user name and password if required.

02. Select Standard Test (when using a site plan) or Quick Test for general sampling. Under standard testing, select the correct site to be tested.

03. Sample your site by “drawing” a 4” x 4” square. With a vertical and horizontal back-and-forth motion “fill” in the square. Consistency is critical.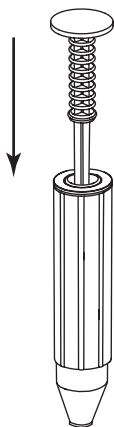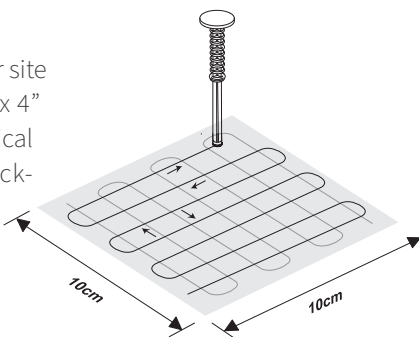


04. Hold the sampler vertically and fully depress to activate.



NOTE: the sampler must be held vertically.

05. Mix gently for two seconds.

NOTE: do not shake.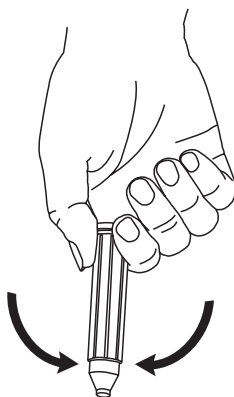


06. With the instrument held upright, press the eject button and insert the sampler.

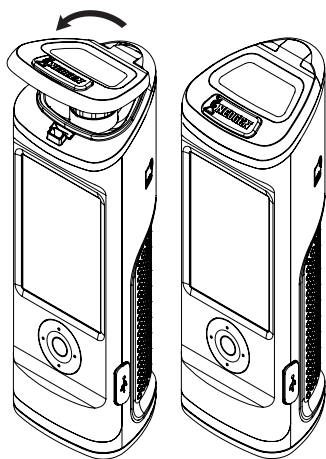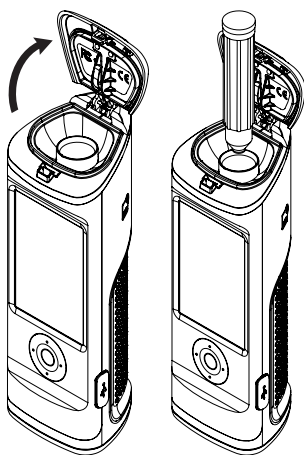07. Depress the sampler all the way down and close the door.
NOTE: keep instrument upright while the sample is being read.

08. The RLU score will display along with a symbol for pass, marginal or fail.

09. On the results screen, users have the option to enter notes for that result, retest that site, or move to the next site in their plan.

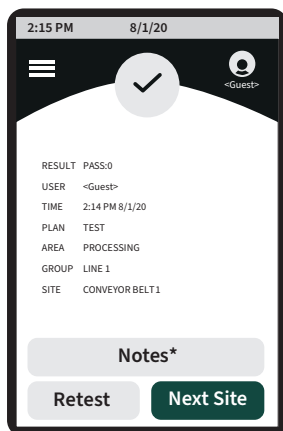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