






## Serial Number

The serial number of the device can be found on a sticker on the left side of the microscope when in front of you (eg. 20200313000 on the illustration below)

### iLine F Pro



## Symbols

---	Direct Current
CE	CE Marking
FCC	FCC Mark, stating the equipment is compliant with the regulations of the US Federal Communications Commission regarding electromagnetic interference
	Do not dispose. The equipment must be properly recycled.
	Safety warning. Read the user manual.
	Power symbol printed on the main instrument switch. Use to turn the instrument on or off.
	Protective conductor terminal
	Earth terminal

## Compliance

The iLine F Pro meets the requirements of the following Directives of the European Union:

- Low Voltage Directive (LVD) 2014/35/EU
- Electromagnetic Compatibility (EMC) Directive 2014/30/EU
- FCC ID: 2AXQY-I3TFMICG02A. *This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.*

## Secure FTP Copy

The remote copy functionality allows you to copy all images captured during your session to a SFTP remote folder.

Enter the SFTP URL and your SFTP credentials. Then, click on *Validate* to ensure that the SFTP information are valid.

See above the “Remote Copy” section for a description of the data copied during the SFTP remote copy process.

## Automatic Clean up

Delete all files that have been safely copied to the remote location. Moreover, files are only deleted if they are older than the specified “delay before cleaning up”.

## Information

This screen displays hardware and software information about the iLine F Pro.

Scroll down to see its serial number and the OsOne software version.

The Use Custom Pump Profile checkbox is for maintenance purposes only.

Camera

Name

Point Grey Research - BFS-U3

Connection

Ready

Serial number

19194745

Status

On

Light Source

Name

N/A - N/A

Connection

Ready

Status

Standby

Wavelength

630nm

Partial coherence

On

Objective Actuator

Name

N/A - N/A

Connection

Ready

Done

Pump

Name

N/A - iLine-F D3HM G02-A Rev. 2.1.0

Controller

Ready

Status

Backup pumping

Use custom pump profile

☐

Cartridge

QR code

I3TF-BCD-S02-B 1.0.5 20190404089

Temperature sensor

Name

N/A - N/A

Connection

Ready

Status

On

About

Done

About

Manufacturer	Ovizio Imaging Systems
Mode	Transmission - D3HM
Model	iLine-F D3HM G02-A Rev. 2.1.0
Serial number	I3TFMIC20200124001
Hardware Version	2.1.0
Firmware Version	1.1.1
OsOne version	7.2.0.22497
Computer Name	I3TF20200124001
IP Address	192.168.178.43

Done

Mode	Transmission - D3HM
Model	iLine-F D3HM G02-A Rev. 2.1.1
Serial number	I3TFMIC20200313000
Hardware Version	2.1.1
Firmware Version	1.1.4
OsOne version	7.3.0.0
Computer Name	ComputerName
IP Address	192.168.1.2

Regulatory

FCC ID	2AXQY-I3TFMICG02A
--------	-------------------

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Done